

Better Pumps for Better Yield!



No Seals, No Bearings, No Particle Contamination!

BPS-200

2.6 bar (37.7 psi)

21 liters/min (5.5 gallons/min)



Pump Casing Levitated Impeller Outlet Rotor-Magnet Motor/Bearing Stator

Figure 1: Schematic of the main elements of the maglev centrifugal pump

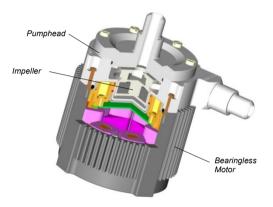


Figure 2: Maglev pump motor with pump head

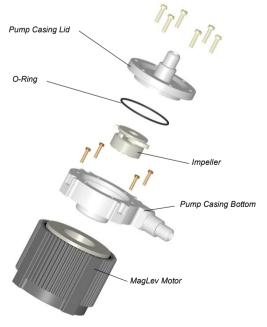


Figure 3: Disassembled pump head

REVOLUTIONARY MAGNETICALLY LEVITATED CENTRIFUGAL PUMP

The *BPS-200* pump system is a revolutionary centrifugal pump that has no bearings to wear out or seals to break down and fail. Based on the principles of magnetic levitation, the pump's impeller is suspended, contact-free, inside a sealed casing and is driven by the magnetic field of the motor (*Figure 1*).

The impeller and casing are both fabricated from chemical-resistant high purity fluorocarbon resins. Together with the rotor magnet they make up the pump head.

Fluid flow rate and pressure are precisely controlled by electronically regulating the impeller speed without pulsation. Closed loop flow or pressure control is easy configurable with the usage of an additional pressure or flow sensor. Various system configurations are possible as explained in the following sections.

SYSTEM BENEFITS

- Extremely low particle generation due to the absence of mechanically contacting parts.
- Increased equipment uptime.
- Lower maintenance costs by eliminating valves, bearings, rotating seals and costly rebuilds.
- Reduced risk of contamination due to the self-contained design with magnetic bearings.
- Very gentle to sensitive fluids due to low-shear design.
- No narrow gaps and fissures where particles or microorganisms could be become entrapped.
- Smooth, continuous flow without pressure pulsation.
- Electronic speed control.
- Compact design compared to pneumatic and mag-drive pumps.
- Proven technology in medical and semiconductor industry (MTBF > 50 years).

APPLICATIONS

- Semiconductor wet processing
- Flip chip and advanced packaging.
- Solar cell production.
- Flat panel display manufacturing.
- Hard-disk fabrication
- Printer ink handling.
- Pharmaceutical production.



STAND-ALONE SYSTEM CONFIGURATION

The stand-alone configuration of the *BPS-200* pump system consists of a controller with an integrated user panel allowing the operator to set the speed manually (*Figure 7*). The speed is automatically stored in the internal EEPROM of the controller. As an option, the speed can also be set with an analog signal (see specification for *Position 3a* in *Table 2*).

EXTENDED SYSTEM CONFIGURATION

The extended version of the *BPS-200* pump system (*Figure 8*) consists of a controller with an extended PLC interface. The PLC interface allows the speed to be set via an external signal, facilitating precise closed-loop flow or pressure control when either a flow or pressure sensor is integrated into the system. A computer can be connected via a USB interface to allow communication with the *Levitronix® Service Software*. Hence parameterization, firmware updates and failure analysis are possible.

Precise ultrapure flow control systems can be realized with the BPS-200 pump system in combination with LEVIFLOW® flowmeters. Levitronix® provides either turnkey solutions for closed-loop flow control or helps to design your own flow control system. Experience has been gained with fluids such as CMP slurries, surface-conditioning chemicals, plating solutions, ultrapure water and solvents. A block-diagram for a typical point of use flow control system with an additional pre-pressure regulator is shown in Figure 5. The versatility of Levitronix® flow control systems goes far beyond the capabilities of simple flow controllers. In addition to the flow control function, the Levitronix® control firmware comes with several condition monitoring features to monitor the integrity of the fluid circuit. Levitronix® flow control systems can generate alarms for preventive filter exchange, no-flow conditions or line clogging. Dynamic Condition Trending (DCT) enables failure prediction and scheduling of preventive maintenance (Figure 6).

ATEX / IECEx SYSTEM CONFIGURATION

Specific Levitronix® motors are ATEX / IECEx certified and allow therefore installation of motor and pump head within an ATEX Zone 2 area (see Figure 9). An Ex conform solution is needed for the motor cables to leave the Ex area. One option is an Ex certified cable system as listed in Table 4 (see Pos. 8) and shown in Figure 13.

- ATEX / IECEx certified for Category 3G/D (Zone 2 for Gas and Zone 22 Dust).
- Thermal classification T5 (< 100 °C = 212 °F) for maximum liquid temperature of 90 °C / 194 °F.</p>
- ATEX / IECEx marking of motor with pump head:
 - CE II 3G Ex nA IIC T5 Gc
 - C€ E II 3D Ex tc IIIC T100°C Dc

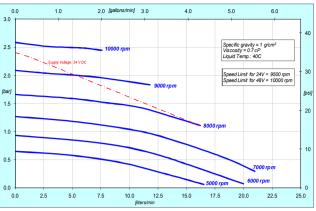


Figure 4: Pressure/flow curves for aqueous liquids (similar to water)

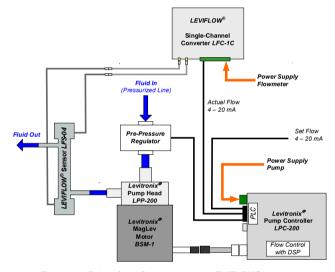


Figure 5: Point-of-use flow control with LEVIFLOW® flowmeter LFS-04 and pump system BPS-200

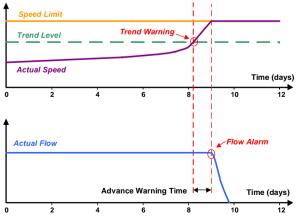


Figure 6: Dynamic Condition Trending (DCT)



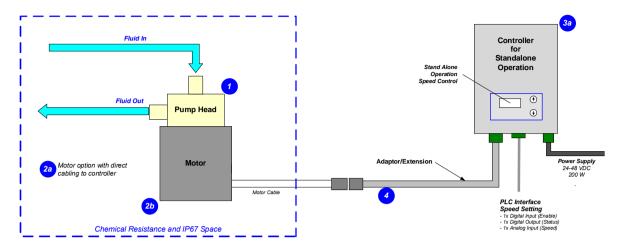


Figure 7: System configuration for standalone operation (Speed setting with integrated user panel)

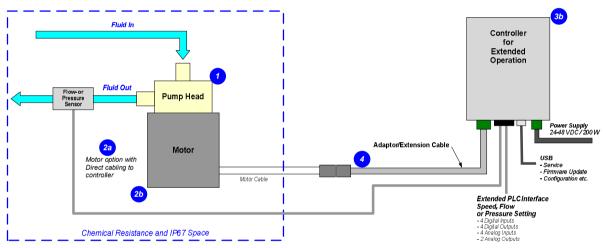


Figure 8: Extended operation (flow or pressure control) with extended controller

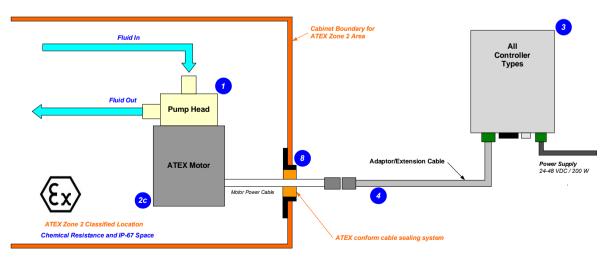


Figure 9: System Configuration for ATEX / IECEx applications



DIMENSIONS OF MAIN COMPONENTS

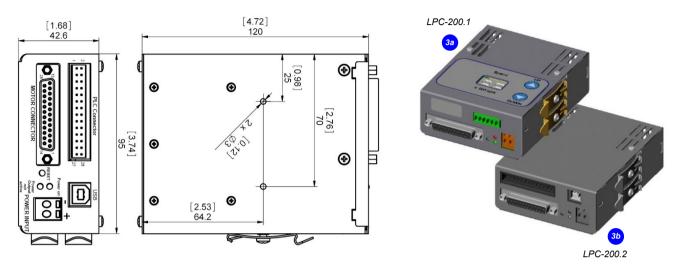


Figure 10: Dimensions of controllers LPC-200.2 (same basic dimensions for LPC-200.1)
(dimensions in mm and [inch])

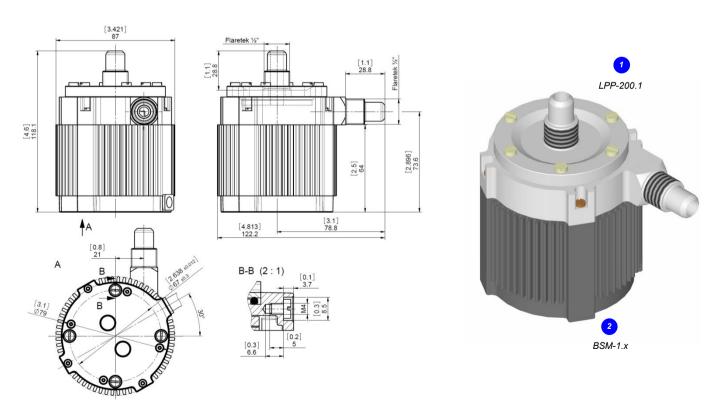


Figure 11: Dimensions of motor with pump head (Dimensions in mm and [inch].)



System Name	Article #	Pump Head	Motor	Controller	Note
BPS-200.3/4	100-901 93 / 94	LPP-200.1	BSM-1.4	LPC-200.1/2	Direct cabling between motor and controller.
BPS-200.1/2	100-901 91 / 92		BSM-1.3	LPC-200.1/2	Adaptor/Extension (0.5 - 10m) cables according to Table 3 have to be ordered as separate article.
BPS-200.14 / 15 (ATEX)	100-905 56 / 57	(High Viscosity and Density)	BSM-1.6 (ATEX)	LPC-200.1/2	Adaptor/Extension (0.5 - 10m) cables according to Table 3 have to be ordered as separate article. ATEX Cable Sealing System can be ordered according to Table 4. ATEX and IECEx certified.
BPS-200.37 / 24	100-907 73 / 03	LPP-200.11	BSM-1.4	LPC-200.1/2	Direct cabling between motor and controller.
BPS-200.40 / 41	100-907 76 / 77	(Improved bubble and	BSM-1.3	LPC-200.1/2	Adaptor/Extension (0.5 - 10m) cables according to Table 3 have to be ordered as separate article.
BPS-200.44 / 45 (ATEX)	100-907 80 / 81	priming behavior)	BSM-1.6 (ATEX)	LPC-200.1/2	Adaptor/Extension (0.5 - 10m) cables according to Table 3 have to be ordered as separate article.
BPS-200.50 / 51 (ATEX)	100-911 10 / 11	LPP-200.10 (Antistatic)	BSM-1.6 (ATEX)	LPC-200.1/2	ATEX Cable Sealing System can be ordered according to Table 4. ATEX and IECEx certified.

Table 1: Standard system configurations
Note 1: All systems are CE, IECEE CB scheme and ETL (NRTL) certified.

Pos.	Component	Article Name	Article #	Characteristics	Value / Feature	
1a		LPP-200.1 (High Viscosity/Density) LPP-200.11 LPP-200.10 (Antistatic) (Improved priming/bubble behavlor)	100-90281 100-90702 100-90684	Impeller / Pump Housing Sealing Ring Fittings	LPP-200.1/11: High Purity PFA / PTFE LPP-200.10: PFA / antistatic PTFE Kalrez® perfluoroelastomer 1 LPP-200.10: Pillar FM ½" LPP-200.10: Pillar FM ½"	
1b 1c	Pump Head			Max. Flow Max. DiffPressure Max. Viscosity / Density Max. Liquid Temp	21 liters/min / 55 gallons/min 2.6 bar / 37.7 psi LPP-200.11/10: 50cP / 1.8 g/cm3 LPP-200.11/10: < 20 cP 90°C / 194°F	
				Pumphead Selection	LPP-200.1: Usage for higher viscosity and density LPP-200.11: Usage for aqueious liquids and improved priming/bubble robustness	
2a	Motor	BSM-1.4	100-10005	Housing	ETFE (chemical resistant) coated Aluminum, waterproofed (IP67)	
Za				Cable / Connectors	1x 5m cables with FEP jacket / 1x D-SUB (direct cable to controller)	
2b	Motor	BSM-1.3	100-10004	Cable / Connectors	1x 2m cable with FEP jacket / 1x circular (M23, IP-67) (needs extension adaptor cable for connection to controller)	
2c	Motor (ATEX / IECEx)	BSM-1.6	100-10063	Cable / Connectors	1x 2m cable with FEP jacket / 1x circular (M23, IP-67) (needs extension adaptor cable for connection to controller)	
	motor (/ 11271/ 1202X)			ATEX / IECEx Marking	(€ 1 3 G Ex nA IIC T5 Gc	
3a		LPC-200.1	100-30008 (Controller with Enable connector incl. in 100-90335)	Electrical Power / Voltage	200 W / 24 - 48V DC	
	Standalone Controller (User Panel)			Interfaces for Standalone Controller	Panel to set speed (automatic storage on internal EEPROM)	
					1x analog input ("Speed") 4 - 20 mA PLC with 1x digital input ("Enable") 0 - 24 V (optocoupler) 1x digital output ("Status") 0 - 24 V (relais)	
				Standard Firmware	C2.25	
3b	Extended Controller (PLC and USB)	LPC-200.2	100-30009 (Controller with PLC connector incl. in 100-90336)	Interfaces for Extended Controller	- up to 4 digital inputs 0 - 24V (optocoupler) - up to 4 digital outputs 0 - 24V (relais) PLC with up to 2 analog inputs 4 - 20mA - up to 2 analog inputs 0 - 10 V - up to 2 analog outputs 0 - 5 V	
	(i LO and OOD)				USB interface (for service and system monitoring)	
				Standard Firmware	C2.48	

 Table 2: Specification of standard components

 Note 1: Kalrez® is a registered trademark of DuPont Dow Elastomers
 Note 2: Pump heads with Pillar FM ½" fittings available.

Pos.	Component	Article Name	Article #	Characteristics	Value / Feature
4	Extension Adaptor Cable for Power	MCA-1.4-05 (0.5m) MCA-1.4-30 (3m) MCA-1.4-50 (5m) MCA-1.4-70 (7m) MCA-1.4-100 (10m)	190-10228 190-10170 190-10187 190-10229 190-10230	Jacket Material Connectors	PVC-jacket Circular Hummel to D-SUB connector

Table 3: Specification of adaptor/extension cables

Pos.	Component	Article Name	Article #	Characteristics	Value / Feature	
5	Air Cooling Module	ACM-1.1	190-10003	Material / Connection Port Air Pressure / Consumption	PVDF / NPT '/." ~ 0.5bar (7.2 psi)	
6	Mounting Base Plate	MBP-1.1	190-10004	Material	PVDF	
7 (a,b,c, d,e)	Impeller Exchange Kits	IEK-200.1 (for LPP-200.1) IEK-200.2 (for LPP-200.11) IEK-200.8 (for LPP-200.10)	100-90517 100-90524 100-91109	Impeller (a) / Sealing O-Ring (b) Pump Casing Screws (c) Pump Motor Screws (d) Exchange Tool IET-1.1 (e)	PFA / O-Ring, Kalrez, 50.52 x 1.78 IEK-200.1/2: 6 pcs M4 x 15 mm, PVDF IEK-200.8: 6 pcs M4 x 16 mm, PTFE coated stainless steel IEK-200.8: 4 pcs M3 x 16 mm, PVDF IEK-200.8: 4 pcs M3 x 14 mm, stainless steel (SS) POM-C	
8	ATEX Cable Sealing System	ACS-A.1	100-90292	Sleeve (a) / Gasket (b) Frame (c) Cable Module (d)	Stainless Steel and EPDM Roxylon (EPDM rubber) Roxylon (EPDM rubber)	Note: Lubricant (e) and measurement plates (f) are included.
9a	AC/DC Power Supply	TSP 360-124 (Standard Traco Supply)	100-40016	Voltage / Power Output / Voltage Input Dimensions	24 VDC / 360 W 85 – 132 / 187 – 264 VAC, autoselect 125 x 125 x 80 mm	
				Certification	UL, CSA, CB, Semi F47	
9b	AC/DC Power Supply	TSP 360-148	100-40017	Voltage / Power Output	48 VDC / 360 W (other specifications same as 9a)	
10	Screw Set	Screw Set SS+PTFE	100-90447	Number/Dimensions Material	4 pcs M3x14 and 6 pcs M4x16 Stainless Steel+ PTFE coating	Needed to increase maximum static pressure of pumphead.

Table 4: Specification of accessories













Figure 12: Pump system BPS-200 with main standard components











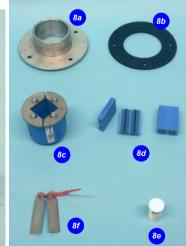


Figure 13: Accessories



LEVITRONIX® THE COMPANY

Levitronix® is the world-wide leader in magnetically levitated bearingless motor technology. Levitronix® was the first company to introduce bearingless motor technology to the Semiconductor, Medical and Life Science markets. The company is ISO 9001 certified. Production and quality control facilities are located in Switzerland. In addition, Levitronix® is committed to bring other highly innovative products like the LEVIFLOW® flowmeter series to the market.



Headquarter and European Contact

Levitronix GmbH Technoparkstr. 1 CH-8005 Zurich Switzerland

Phone: +41 44 445 19 13
Fax: +41 44 445 19 14
E-Mail: salesEurope@levitronix.com

US Contact

Levitronix Technologies Inc. 10 Speen Street, Suite 102 Framingham, Massachusetts 01701 USA

Phone: +1 508 861 3800 Fax: +1 508 861 3837 E-Mail: salesUS@levitronix.

Japan Contact

Levitronix Japan K.K. Wing Eight 5floor, 4-16-4 Asakusabashi, Taito-ku Tokyo, 111-0053 Japan

Phone: +81 3 5823 4193 Fax: +81 3 5823 4195 E-Mail: salesJapan@levitronix.com

Taiwan Contact

Levitronix Taiwan 5F, No. 251, Dong Sec. 1, Guangming 6th Rd., Chu Pei City, Hsin-Chu 302, Taiwan, R.O.C.

Phone: +886 3 657 6209
Fax: +886 988 321472
E-Mail: salesAsia@levitronix.com

PL-4009-00, Rev07, DCO# 20-144

First Release: 03-Oct-2008

Last Update: 09-Jul-2020